

GETTING STARTED WITH OXYGEN OPTODES (OP-SERIES)

1. UNPACKING

- Remove the grey shock-absorbing plastic net and inspect the sensor visually. Leave the sensor in the protection tube for testing.

2. CONNECT THE OPTODE TO THE MICROOPTODE METER

- Remove the cap from the MicroOptode connector
- Remove the cap from the MicroOptode Meter
- Insert the MicroOptode cable into the MicroOptode meter by gently turning clockwise until the plug is locked firmly.

3. CALIBRATE THE SENSOR

- Use air saturated water as one calibration point. This is easily done using the CAL300 calibration chamber. The sensor may be dipped directly into the calibration chamber or the air saturated water may be injected into the protection tube using the calibration cap (see picture).
- Use the Unisense zero O₂ solution as the second calibration point. Inject the zero O₂ solution into the calibration cap and wait for the sensor to respond.
- For alternative calibration method, see the Oxygen MicroOptode Manual.

4. APPROVE THE SENSOR

- Compare the calibration values to Unisense Standard specifications (incl. in sensor box). If necessary, see Troubleshooting in the optode manual or contact support (see below)

5. ADJUSTING THE MEASUREMENT FREQUENCY

- The lifespan of the optodes is determined by the amount of measurements. In order to prolong the lifespan of the optode, adjust the measuring frequency as required for your application
- The frequency can be adjusted in the SensorTrace software

6. STORAGE


- When not in use unplug the optode and store it with the protection tube mounted at room temperature (preferable in low light or darkness).



Unisense
MicroOptode



CAL300 with
microsensors and
bubbling with air.

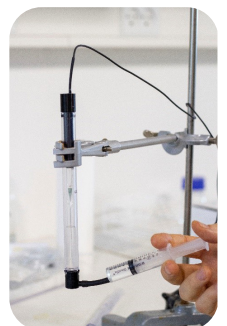

Injecting calibration
liquid into protection tube
using the calibration cap.

USEFUL TOOLS



For support go to
www.unisense.com/support/ or
contact sales@unisense.com



Get the full manuals for all
sensors, equipment & software at
www.unisense.com/manuals/.



O₂ Microsensor
Manual



Calkit-O2
Manual



SensorTrace Suite
Manual



Find SDS for
Calibration Kit here