

Instructions for correct packing of microoptodes

1. Placing the first foam inlay

Place two pieces of foam inlay in the box.



2. Place optode on foam

Place the optode(s) with protection tube(s) and the entire cable on the foam inlay.

Make sure the cable is not bent in any way, and that nothing presses down on it, as this will damage the fiber inside.

The minimum diameter for the rolled up cable is 10 cm.



3. Sensor between the foam inlays

Place the last two pieces of foam inlays on top of the optode(s) and cable(s) to fixate them in the box.



4. Close and ship

Close the lid and push the lid tabs into the slots in the box sides.

